

Product Specification

Product : Antibody Enhancing Solution
Product Number : BDANS250
Package : 250 mL
Storage : RT

PROTOCOL

1. Blocking (with Fast Blocking Buffer).
 2. Washing with PBST or ddH₂O 2 times .
 3. Adding 1st Ab in Antibody Enhancing Solution dilution and incubating at RT for 1~2hr.
 4. This product can effectively enhance the interaction between 1ST Ab and antigen on membrane.
 5. Washing with PBST or TBST 3 times.
 5. Add 2ndAb in lab home buffer for RT 1hr.
 6. Washing with PBST/TBS 3times.
- p.s. The 1stAb in Antibody Enhancing Solution can be reused for 3 times or more.

STORAGE

RT

